



Laboratory Equipment Manufacturer
www.mrclab.com



INSTRUCTION MANUAL FOR **HOP- 1212** LABORATORY HOTPLATES



PLEASE READ THIS MANUAL CAREFULLY BEFORE OPERATION

3, Hagavish st. Israel 58817 Tel: 972 3 5595252, Fax: 972 3 5594529 mrc@mrclab.com

MRC.VER.02-11.10

INSTRUCTIONS for use of LABORATORY HOTPLATES

Model HOP- 1212

230V/ 50Hz, 500W

HEATING SYSTEM :

The Hotplate is actuated by the Temperature Control Knob on the front of the instrument, which regulates heat of the top plate from ambient temperature up to 300⁰C. The pilot lamp indicates the thermostat's operation.

WARNING :

- * the plate may be hot to touch, even when the pilot lamp is off
- * the plate can be kept clean by covering it with Aluminium foil.
If so, cover plate only, and not radiation shields underneath the plate.
- * the instrument should not be used where danger of explosion exists.